

DLAB




**Real-Time PCR System
Accurate 96**



Real-Time PCR System




DLAB Accurate 96 is based on global vision of product design concepts and manufacturing processes. It creatively combines Fresnel lens optical signal acquisition technology, time-resolved signal separation technology and unique temperature control technology. And it reaches international advanced level in sensitivity, multi-color crosstalk, temperature uniformity and accuracy. It supports the application of all common QPCR detection modes.

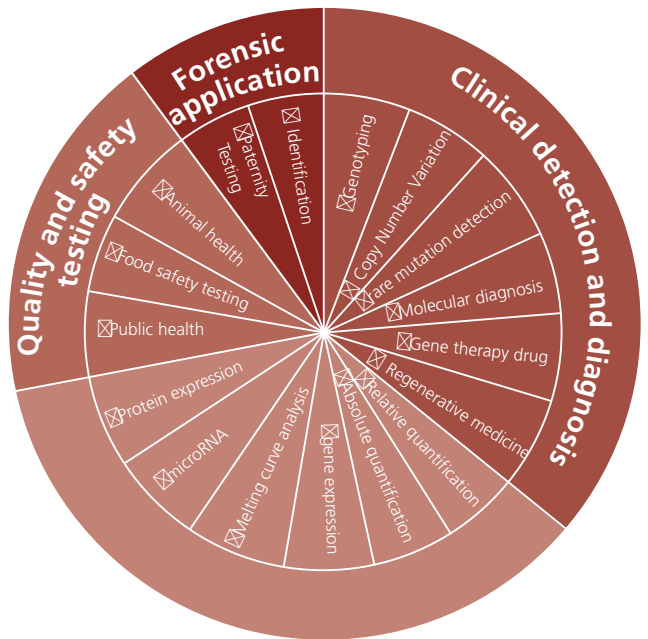

Up to 6 fluorescence detection channels allowing multiplex PCR — Simultaneous detection of 5 target genes in 96 samples

| Channel 6 | Channel 5 | Channel 4 |
|--------------------------|-------------------|------------|
| NED™ /Cy3/ TAMRA™ | ROX™ /Texas Red® | FAM™ /SYBR |
| VIC® /HEX™ /TET/ JOE™ | Cy5® /Quasar® 670 | FAM™ /SYBR |
| Channel 3 | Channel 2 | Channel 1 |

FAM,JOE,VIC,HEX,NED,TAMRA and ROX are trademarks of Life Technologies Corporation and its subsidiaries. VIC and Texas Red are registered trademarks of Life Technologies Corporation and its subsidiaries. Cy5 is a registered trademark of GE Healthcare Bio-Sciences Corp. Quasar are registered trademarks of Biosearch Technologies, Inc.

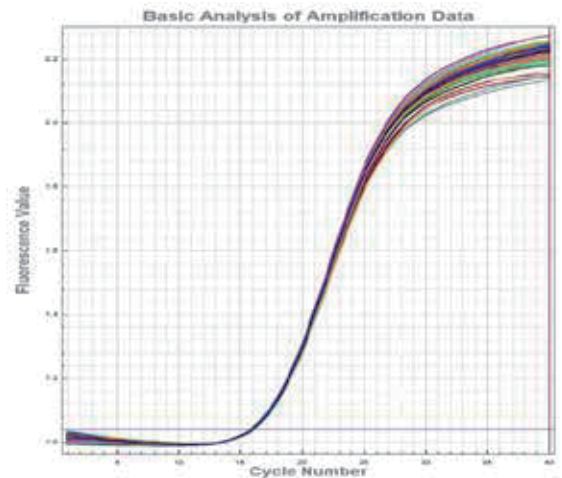



Simultaneous scanning of the six-channel shows that the standard deviation of the Ct value of the FAM channel is <0.07. No fluorescence signal in other channels.

Technical Innovation 1

Effectively reduce multi-color crosstalk and edge effect, no ROX correction required


The multi-color crosstalk caused by the small sample spacing of 96 or 384-well plates has a great influence on the accuracy of the experimental results, especially in multiplex qPCR detection.

The new optical signal detection system and unique time-resolved scanning can reduce non-target sample optical signal collection. Thereby high repeatability of single fluorescent channel can be ensured.

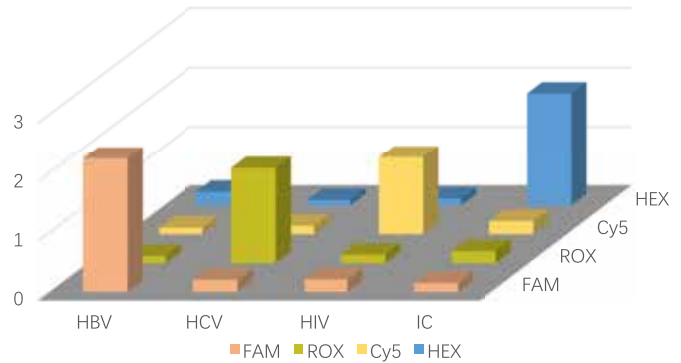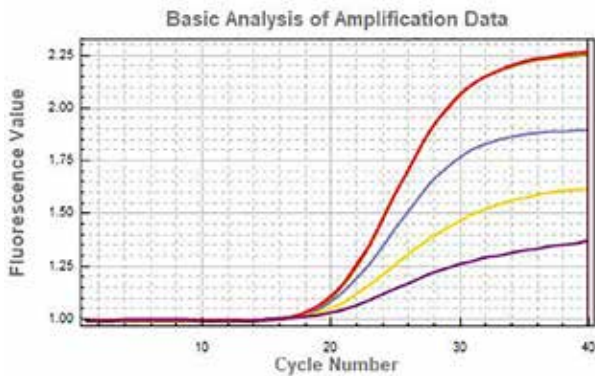


Four different target genes (2 repeats) of FAM/HEX/ROX/Cy5 were simultaneously detected in one reaction tube, and the results showed that there was almost no cross-interference between the different channels.

New optical scanning detection system — High sensitivity /resolution

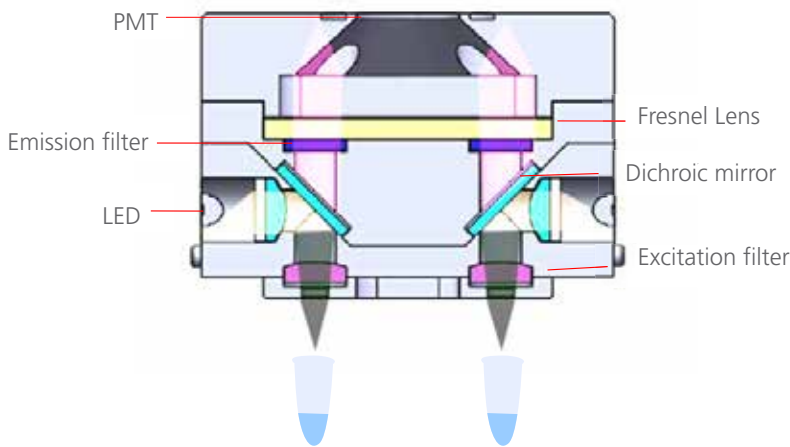

Sectional view of optical detection

- LED light source Efficient and maintenance-free
- Fresnel Lens It greatly reduces the light collection of the non-target area. And the relative position of the detector to the block hole ensures that one optical detection channel is aligned with one target to be tested at the bottom.

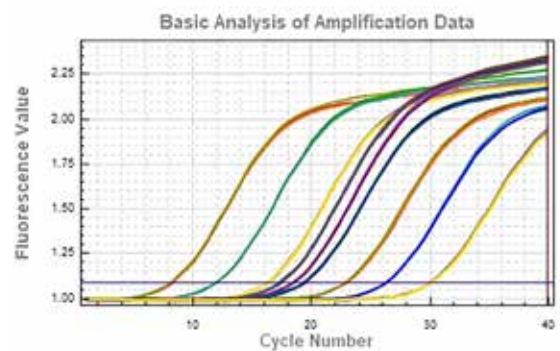


Different concentrations of plasmids were amplified by probe assay (concentration from left to right is 5 μ g, 500 ng, 50 ng, 20 ng, 10 ng, 5 ng, 500 pg, 50 pg, 5 pg), three replicates for per concentration. The Ct values difference of the 10-fold dilution is exactly 3.3. The Ct values difference of the 2-fold dilution is exactly 1.

Real-Time PCR System


Innovative scanning method and time-resolved signal separation technology—High accuracy

- Unique time-resolved scanning method
 - ☒ The different fluorescence signals of the same sample are collected at different times.
 - ☒ The high-speed stepping motor and the highly sensitive detector ensure that all signal acquisition of the entire sample plate is completed in a short time.

- Innovative detection channel arrangement Interlaced arrangement of upper and lower channels further reduces inter-hole and multi-color fluorescence crosstalk.
- Double FAM scanning for melting curve Scan time is shortened a lot.

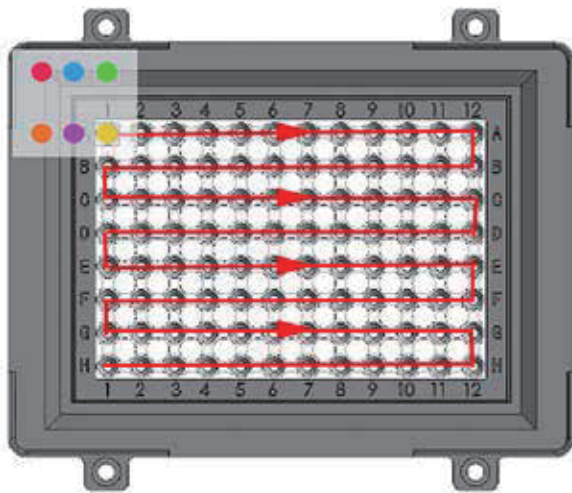


Multi-channel scanning for probe assay

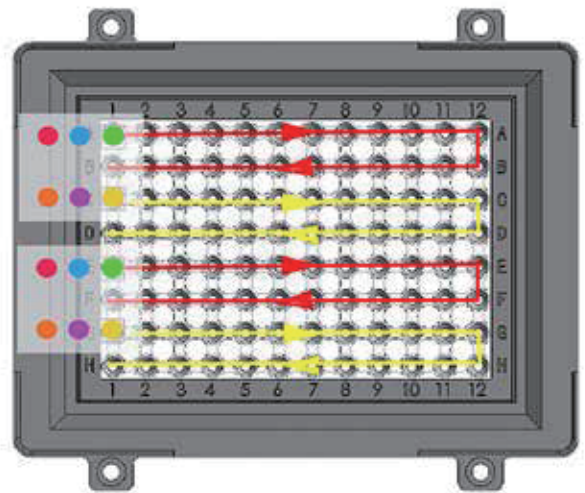

Double FAM scanning for melting curve

Technical Innovation 2

—Unique edge temperature compensation technology

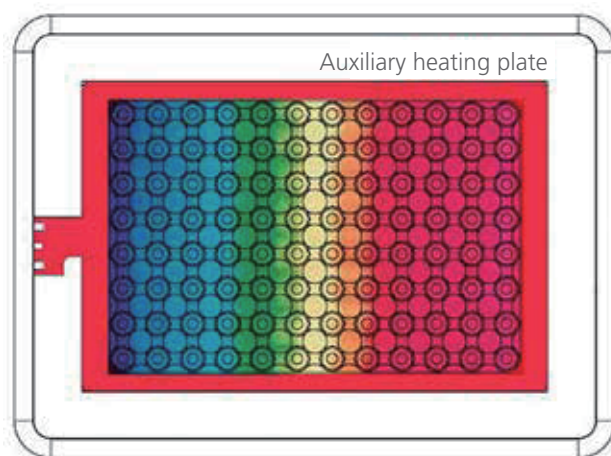

- Both temperature accuracy and uniformity are $\pm 0.2^{\circ}\text{C}$
- Module maximum ramp rate is $6^{\circ}\text{C} / \text{sec}$
- The average sample ramp rate is $2.2^{\circ}\text{C} / \text{sec}$
- With a unique outframe protection design, the unit can achieve very even temperature across the whole plate. Effectively reduce the edge temperature variance.







The T_m values of 96 replicate wells were detected. The results showed that the temperature uniformity difference was $< \pm 0.2^\circ\text{C}@81^\circ\text{C}$.

Temperature profile of 96-well plate



Temperature ($^\circ\text{C}$)

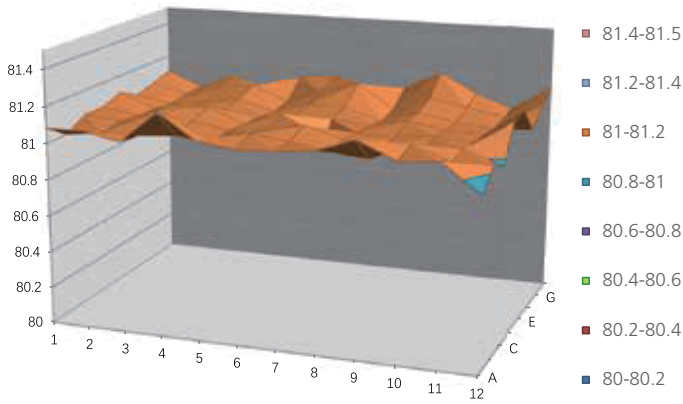

T_m value normal distribution map of 96-well plate

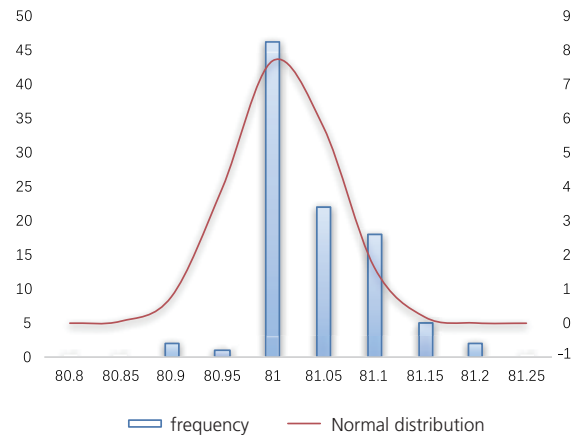

The melting curves of 48 replicate samples

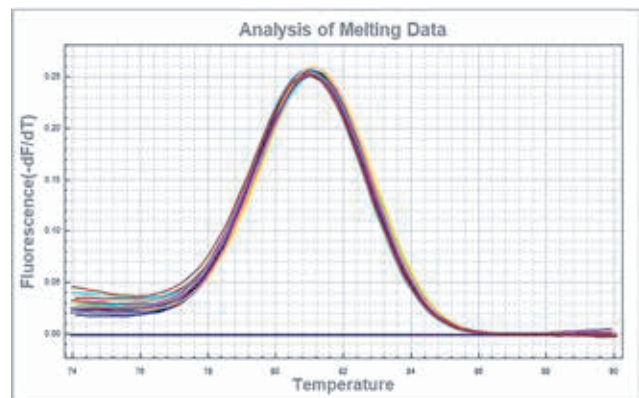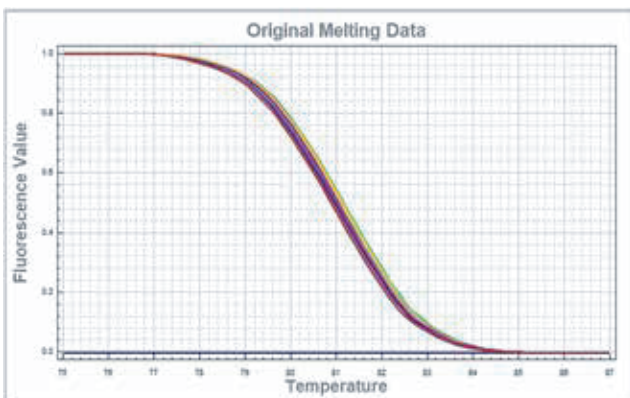


Real-Time PCR System

User-friendly software



Free software upgrades

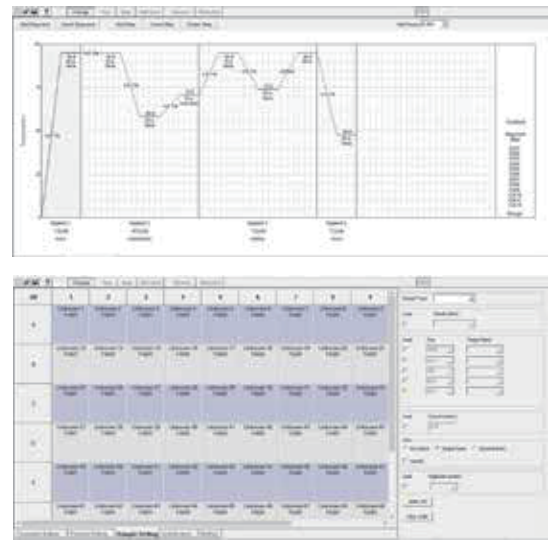

The software can provide you with a full range of solutions for sample detection and result analysis.


One-stop process and solution




Sample preparation and nucleic acid extraction



Prepare PCR mixture

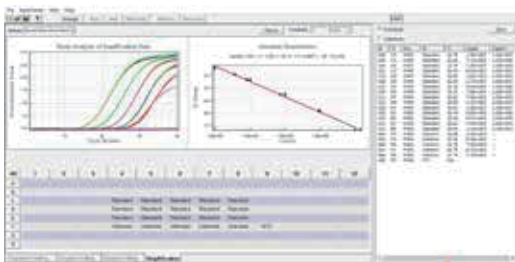

Data analysis

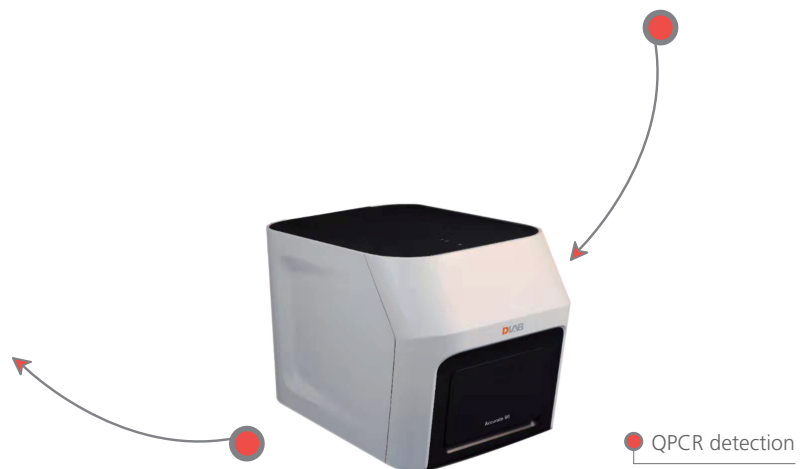

Classic Examples

— Swine fever virus ASFV detection

QPCR was used to detect the swine fever virus DNA in vitro for clinical diagnosis of suspected diseased pigs.

Kit: Swine fever virus (ASFV) nucleic acid detection kit (PCR probe assay) 、 DNA extraction kit (Spin column method)

Method: 5 mL of blood was extract from a live pig syringe to be examined. DNA was extracted by DNA extraction kit. According to the ASFV detection kit operation method, the extracted DNA, the positive control substance, and the negative control substance is separately added to the PCR reaction solution and the enzyme. Then the mixture is centrifuged and tested on real-time qPCR. The reporter and quencher of the TaqMan probe is FAM and TAMRA.


Judgment basis: Positive: $Ct \leq 35$, and the amplification curve has a significant exponential growth curve. Negative: $Ct > 37$, or no significant amplification of the curve. Recommended retest: $35 < Ct \leq 37$.

Data analysis: The Ct value of the positive control is 20, and the negative control product has no obvious amplification curve. The Ct values of the sample 1 and 2 to be tested is 26 and 30 respectively. According to the judgment basis of the kit, the samples to be tested are ASFV positive.

Program:


| Step | Temperature | Time | Cycle | End point |
|------|-------------|-------|-------|-----------|
| 1 | 95°C | 10min | 1 | No |
| 2 | 94°C | 15sec | 40 | No |
| | 55°C | 30sec | | Yes |

Amplification curve:

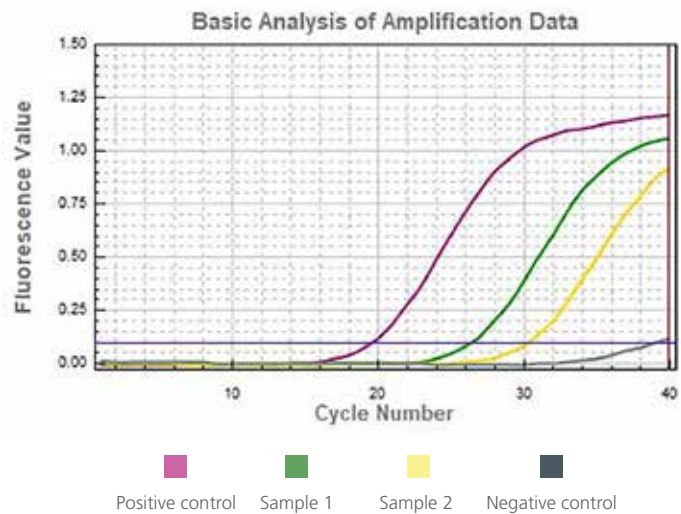

Technical Parameters

| Temperature control system | | Detection system | |
|---|---|---------------------------------------|--|
| Sample capacity | 0.1ml PCR tubes×96, 8×12 PCR plate or 96 well plate × 1 | Excitation light source | 4/6 monochrome high efficiency LEDs |
| Reaction volume | 10-50 μl | Detection device | PMT |
| Thermal cycle technology | Peltier | Detection mode | Time-resolved signal separating technology |
| Max. Heating/Cooling rate | 6.0° C/s | Excitation/detection wavelength range | 455-650nm/510-715nm |
| Heating temperature range | 4 – 100 °C | Fluorescent channels | 4/6 channels |
| Temperature accuracy | ± 0.2°C | Supported dye | FAM/SYBR Green, VIC/JOE/HEX/TET, ABY/NED/TAMRA/Cy3, JUN, ROX/ Texas Red, Mustang Purple, Cy5/LIZ |
| Temperature uniformity | ±0.2°C @60°C , ±0.2°C @95°C | Sensitivity | Single copy gene |
| Temperature gradient setting range | 30–100°C | Resolution | 1.33 folds copy number difference can be distinguished in single-plex qPCR |
| Temperature gradient difference setting range | 1 – 36°C | Dynamic range | 10 orders of magnitude copies |

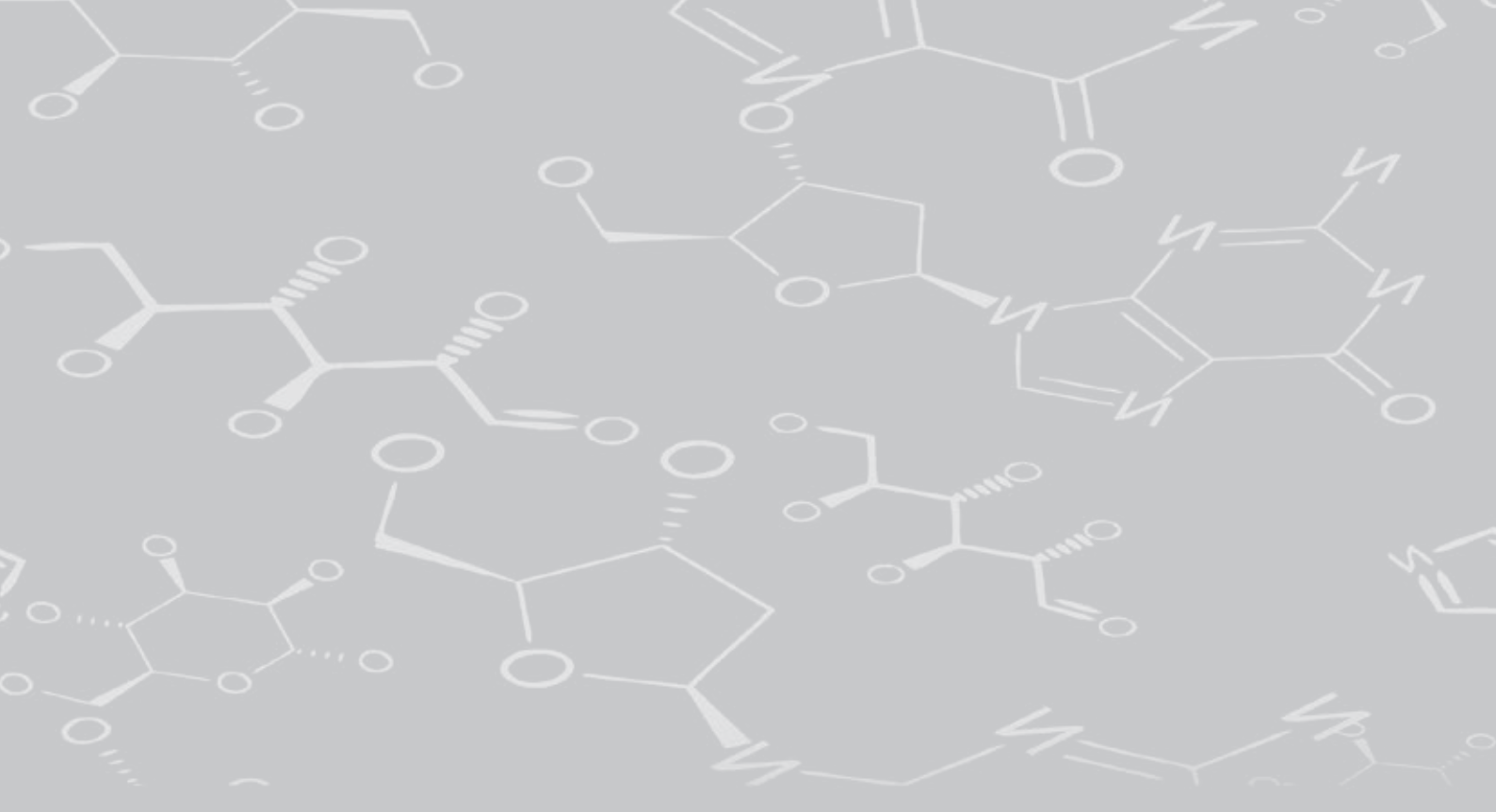

DLAB Scientific Inc.

Add: 775 Rivera St, Riverside, CA 92501, USA

Tel: +1- 747- 230-5179

Fax: +1-909-230-5275

E-mail: info@dlabsci.com

Service contact: service@dlabsci.com

web: www.dlabsci.com

