



## MicroPad Staphylococcus aureus Count Plate

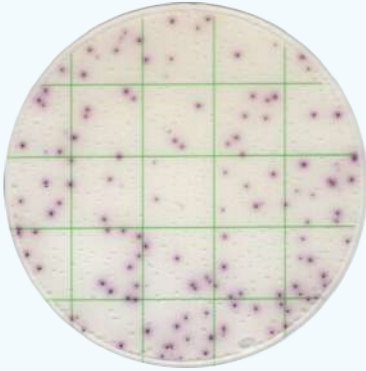


**Product Code: MCPG00225**    **25 plates / pack**

Staphylococcus aureus Count Plate is a sample-ready-culture medium system which contains readymade Baird-Parker medium, a cold-water-soluble gelling agent indicator and selective inhibitor. It can be used in direct counting of staphylococcus aureus colonies after 24h incubation. The result is consistent with ISO standards and commercial counting plates.


**Read the user manual carefully before test, and follow the instructions.**  
**Failure to do so may lead to inaccurate results.**

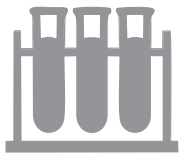


**Sample Preparation**

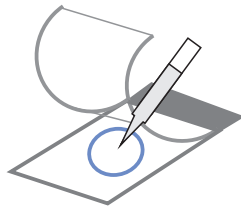

**Sampling**

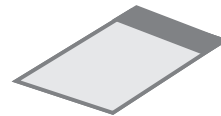

**Incubation**

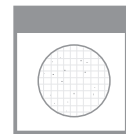

**Enumeration**

Incubate the plates at  $36 \pm 1$  °C for  $24 \pm 2$  hours.

For aquatic products, please incubate at  $30 \pm 1$  °C for  $72 \pm 3$  hours for. Up to 6 plates can be stacked in one incubation holder. On this Staphylococcus aureus count plate, Staphylococcus aureus is red colonies, other staphylococcus is light pink green or colorless colonies. Majority of gram-negative and other gram-positive bacterium cannot grow in this plate, or they can present as blue colonies.